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COMPRESSED AIR THERMODYNAMICS

Compressed air is one of the most important, expensive and versatile commodities used in the industries. The study of the compression phase and its associated transient or not phenomena can only be understood if an efficient data collection system is available. Besides, according to the use, it must have specific characteristics, not only related to pressure, but also as concerns the different contaminants such as bacteria, viruses and dusts in the in-taken air, metal dusts and rust particles in the compressor and the tubes, lubricating oil, water condensate.

All creates a mixture that is the main cause of the faults on the installations and that must, if necessary, be removed. This computerized system is designed so that students can collect and utilize operational data, which in normal circumstances should be difficult to obtain manually.

COMPUTERIZED COMPRESSED AIR PRODUCTION AND TREATMENT TRAINER mod. TTDAC/EV

The system consists in two trainers. The first one, connected to the PC, enables to display the compressed air thermodynamic cycle. The second one, provided with refrigeration cycle, enables the compressed air drying. The unit consists of:

- Industrial components and instruments, which are properly connected and operating, easily accessible and visible
- Two-speed motor unit
- Two-stage compressor with pulley
- Personal computer (supplied upon request) for data acquisition and processing from the system
- Compressed air tank unit
- Refrigeration unit
- 2 tanks for heat exchange
- Instruments and sensors connected to the computer
- Safety devices: E.L.C.B. and standard protections.

The equipment is provided with a personal computer (upon request), the connection interfaces, the proper transducers and sensors.

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TRAINING PROGRAM

- Analysis of the pressure/volume diagram of the 1st and 2nd stage with experimental test
- Calculation of the mean effective pressure of the 1st and the 2nd stage and the power absorbed by the compressor
- Use of the air T/S diagram for the experimental determination of the cycles and heat balances characteristics
- Analysis of the compressor efficiency
- Checking the safety devices
- Analysis of the air thermodynamics with detailed examination of dew point and water content
- Specific examination of the compressed air characteristics with use of different filters and refrigerant circuit
- Analysis of the air contaminants and examination of the possible effects on possible uses
- Air side and freon side heat balance of the cycle; yields.

TECHNICAL SPECIFICATIONS

- Color synoptic diagrams, with 11 signaling lamps, 5 taps, diagrams and switches, mounted on painted, baked and finished steel structures
- Electrical instruments (2 voltmeters, 2 ammeters and 3 wattmeters in class 1.5); 2 pressure gauges range 0-10 and 0-30 bar, respectively, accuracy $\pm 0.6\%$ f.s.
- Motor unit: 2.5-KW three-phase (4-8 poles) with thermal protection
- Air intake chamber with filter, direct reading flow meter, accuracy $\pm 3\%$ f.s. Direct reading hygrometer with transducer and interface to computer, accuracy $\pm 2\%$ f.s.
- Two-stage compressor, 369 cc at the 1st stage and 108 cc at the 2nd stage, respectively; connected to a position encoder and a turbine air flow meter interfaced to the computer
- Air analysis outlet chamber with throttle valve and hygrometer
- Refrigerant circuit: 200-W compressor, forced air condenser, capillary pipe, drying filters, humidity indicator, freon/glycol exchanger
- Compressed air circuit with 8 quick acting couplings for connections to either stage of the compressor, 2 air filters with interchangeable cartridges (2 cartridges with efficiency of the 93% and the 99.9% at 0.1 μ respectively), automatic discharge valve with condensate level indicator, 2 glycol/compressed air exchangers, Plexiglas observation chamber with glycol for heat governor
- Digital thermometers with probes, accuracy $\pm 1\%$ f.s.
- 200-l tank complete with accessories
- Control software
- Optional PC.

OPTIONAL ITEMS

- Personal Computer, minimum configuration: IBM-compatible, 16-MB RAM, VGA and SUPER VGA graphic card, serial interface, parallel interface, Windows 95/98.

POWER SUPPLY

400 V - 50 Hz three-phase and 230 V - 50 Hz single-phase
230 V - 60 Hz three-phase and 110 V - 60 Hz single-phase (upon request)

DIMENSIONS AND NET WEIGHT

mod. TTAC/EV: 130x80x180 cm – 154 kg
mod. TDA/EV: 130x80x180 cm – 113 kg
Tank: 60x60x170 cm – 54 kg

THEORETICAL-EXPERIMENTAL TEXTS

- Theoretical-experimental manual for compressed air thermodynamics

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